

LAARMANN[®]

Innovators in Solids

GREEN COFFEE WITH CM100 CUTTING MILL



Laarmann Group B.V.
Op het Schoor 6
6041 AV Roermond
Netherlands

info@laarmann.eu
www.laarmann.eu
webshop.laarmann.eu
+31 6 23 40 00 33

GREEN COFFEE APPLICATION

Product description:

Green Coffee.



884gram Sample milled in cutting mill CM100 with 4 mm sieve and 2mm sieve.



Sieved with a 2 mm sieve 90% smaller than 2 mm.



Grinding system: CM100 Cutting Mill

Speed: 500 rpm

Accessories: Bottom sieve of 4,00 mm / 2,00 mm

Sample material: Green coffee

Product feed size: Original bean

Grinding time: 3 minutes for 884 grams Result after sieving: homogeneously milled mainly < 2,00mm

Recommendation

In order to grind homogeneously green coffee we recommend the use of the CM100 Cutting Mill under the described parameters. With the 4,00 mm sieve it is really easy to grind 1 kg or larger sample quantities up to several kilos.